


de Drucker (PR)	en Printer (PR)	cs Tiskárna (PR)
da Printer (PR)	el Εκτυπωτής (PR)	en-US Printer (PR)
es Impresora (PR)	fr Imprimante (PR)	it Stampante (PR)
nl Printer (PR)	no Skriver (PR)	pt Impressora (PR)

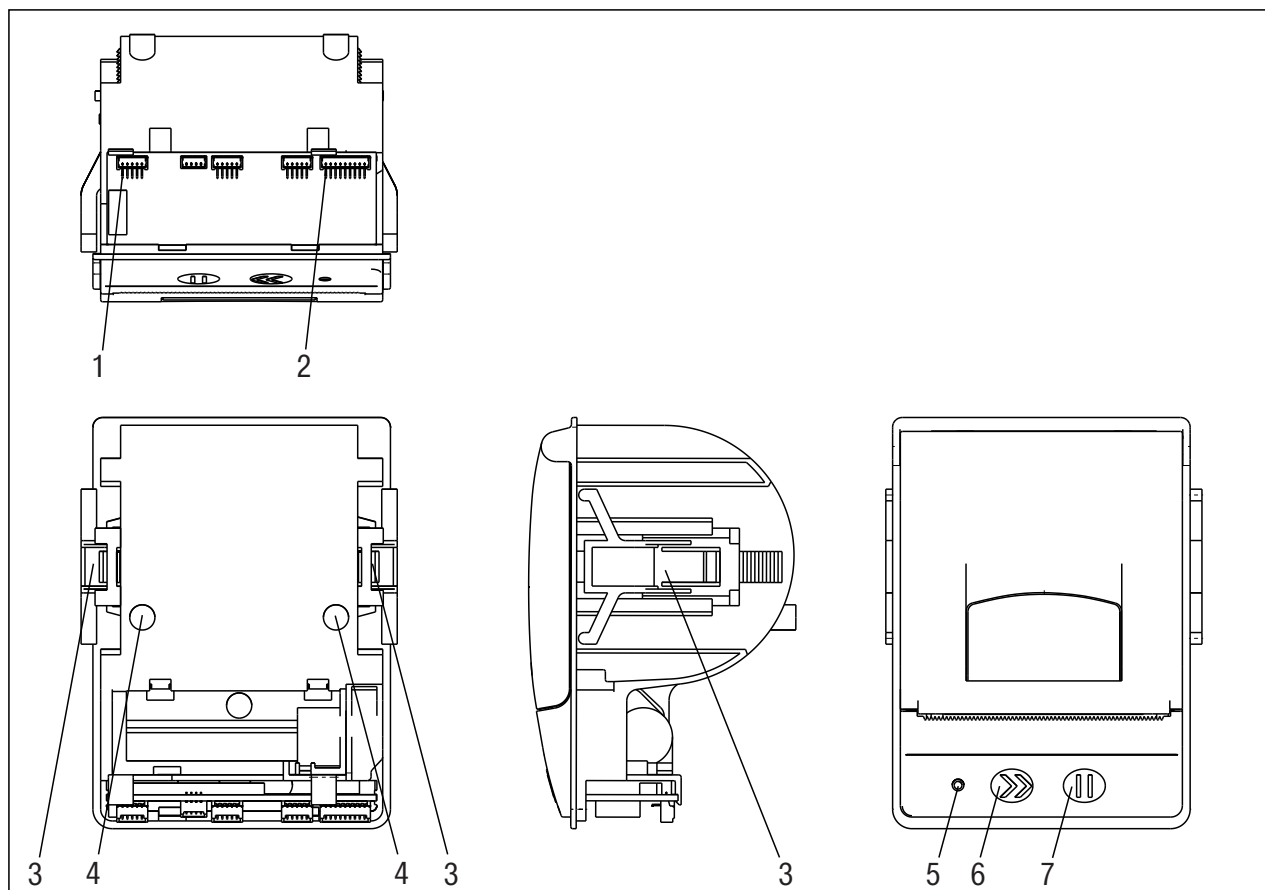

1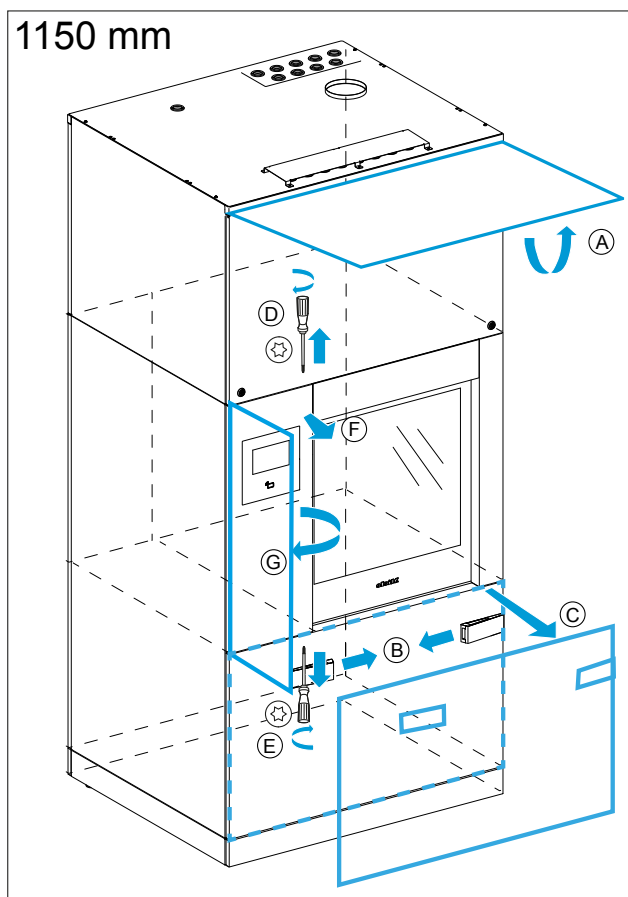


2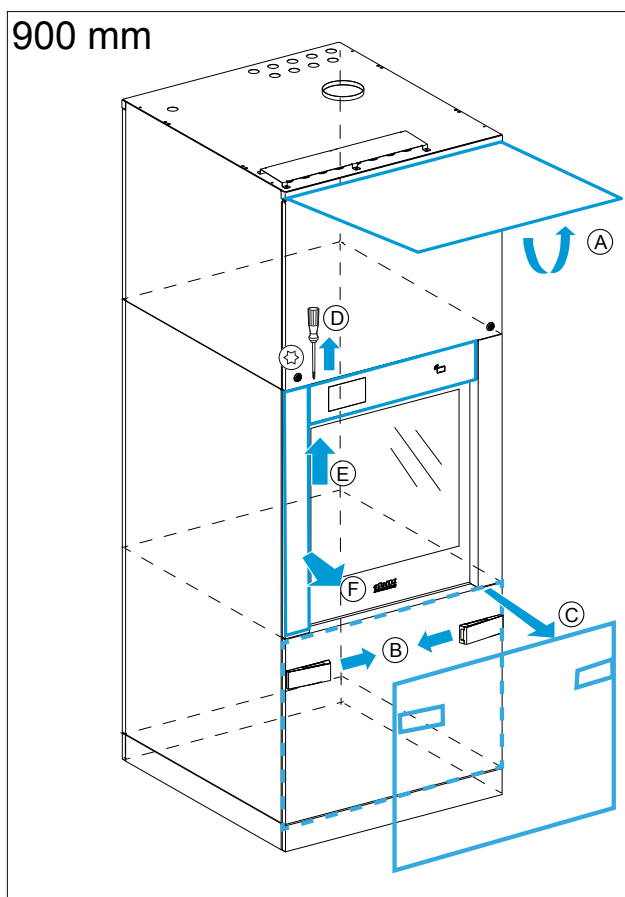


3

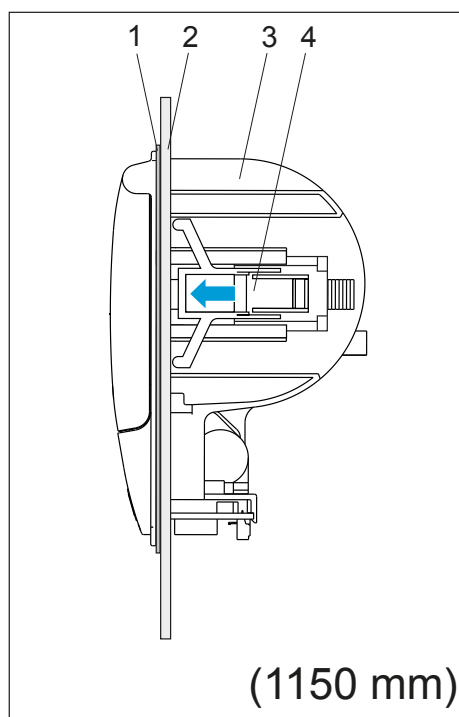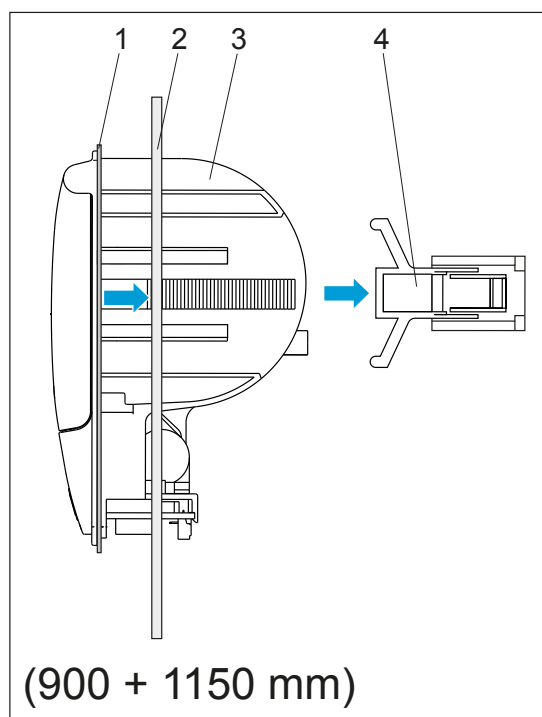
4

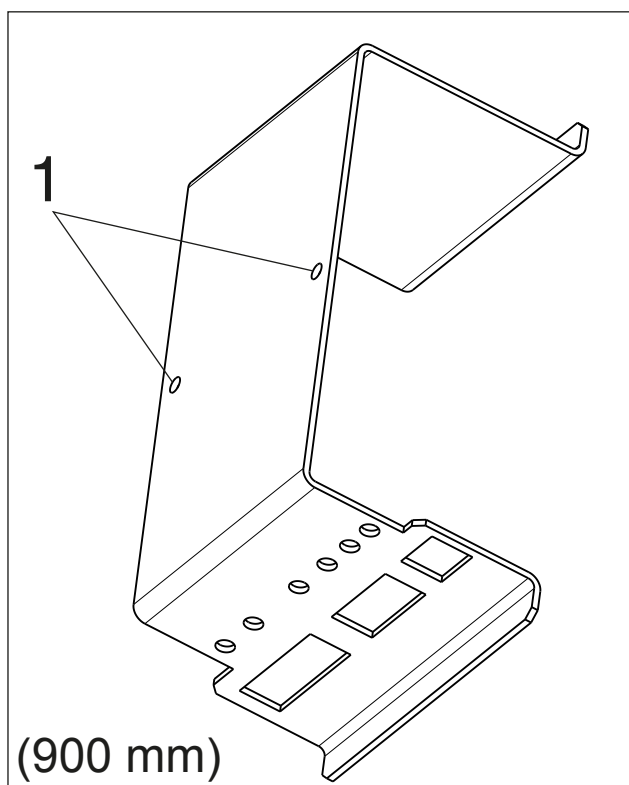5

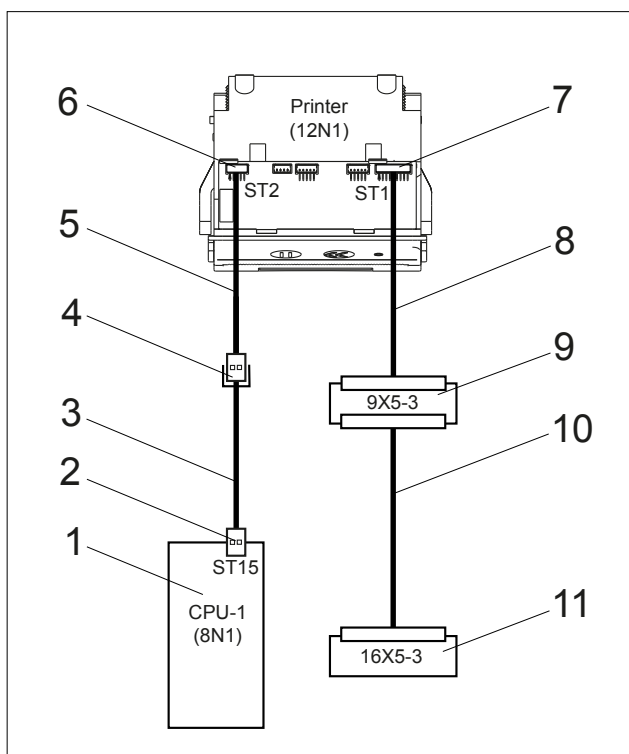6

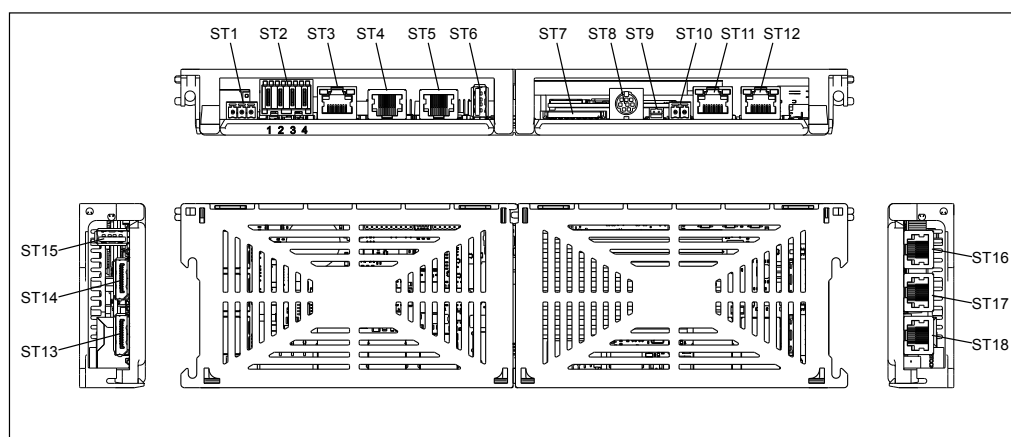7

de

Drucker (PR)

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In order to carry out this work, the appropriate Technical Service Documentation (TSD) is required.

The warning and safety instructions given in the applicable model-specific TSD must be complied with.

Work must be carried out in accordance with the instructions given in the applicable model-specific TSD.

Printer (PR) (12N1)

Legend for Fig. 1:

1. PIN 1 position plug ST2 USB 2.0
2. PIN 1 position plug ST1 power supply 5 V DC / < 3 A
3. Holding clips
4. Thread for fixing screws
5. LED - Operating states
6. Button - Paper feed
7. Button - Confirmation - For continuing with print after paper roll exchange

Printer technology: Thermal, direct

Operational state	Description / State
Constantly off	Power off
1 x flash - Pause ¹⁾	Power on / Not ready
Constantly off	Ready for operation
2 x flash - Pause ¹⁾	Fault: Printer head
3 x flash - Pause ¹⁾	Fault: Paper end or paper feed fault
4 x flash - Pause ¹⁾	Fault: High voltage, low voltage or temperature fault

LED - Operating states

¹⁾ = Pause between the flashing code = 1 s (flashing rhythm = 128 ms on 128 ms off (8 Hz))

Parameters	Unit	Value
Temperature during operation	°C	-5 to 50
Moisture during operation	% uncondensed relative humidity	20 - 85
Storage / Transportation temperature	°C	-40 to 80
Moisture during storage / transport	% uncondensed relative humidity	10 - 90

Ambient parameters

Component	Material
Cover	Sabic Lexan 243R (GE)
Bottom part	Sabic Cycloy C2950

Material**Printer fitting**

Depending on model (900 mm or 1150 mm width) only one of the following work stages can be carried out:

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Subsequently continue with the work step **Printer preparation for fitting**.

Corner piece (wide) opening on unloading side

- ↺ Open the top box panelling service flap (A), see Fig. 2.
- ↺ Release the locks on the bottom part service panel (B), see Fig. 2.
- ↺ Remove the bottom part service panel (C), see Fig. 2.
- ↺ Remove the top fixing screw (D) and the bottom fixing screw (E) from the corner piece (wide), see Fig. 2.
- ↺ Slide out the corner piece (wide) partially from the machine (f), see Fig. 2.
- ↺ Open the corner piece (wide) (G), see Fig. 2.
- ↺ Remove the printer opening cover plate on the corner piece.
- ↺ Continue with the work step **Printer preparation for fitting**.

Corner piece (narrow) removal on unloading side

- ↺ Open the top box panelling service flap (A), see Fig. 3.
- ↺ Release the locks on the bottom part service panel (B), see Fig. 3.
- ↺ Remove the bottom part service panel (C), see Fig. 3.
- ↺ Remove the fixing screw for the corner piece (narrow) and fascia (D), see Fig. 3.
- ↺ Lift the fascia slightly and use a piece of wire or a cable tie inserted through the screw hole to secure it to the top box panelling holder.
- ↺ Lift the corner piece (narrow) slightly (E), see Fig. 3.
- ↺ Remove the corner piece (narrow) (F), see Fig. 3.
- ↺ Remove the printer opening cover plate on the corner piece.

↪ Continue with the work step **Printer preparation for fitting**.

Printer preparation for fitting

↪ Remove the holding clamps (4) from the printer (3), see Fig. 4.

↪ Fit the printer (3) with its seal (1) in the cutout in the corner piece (wide) (2), see Fig. 4.

Fitting printer in corner piece (wide)

↪ Refit the holding clamps (4) in the retainers on the printer (3), see Fig. 4.

↪ Press in the holding clamps (4) until they engage in the corner piece (wide) (2) and thus secure the printer (3) to the corner piece (wide), see Fig. 4.

Fitting printer in corner piece (narrow)

The printer holding clamps are only required if the printer is fitted in the corner piece (wide).

For fitting the printer in the corner piece (narrow) a special holding plate is supplied with the kit.

↪ Fit the holding plate over the printer such that the openings for the connection plug are on the underside.

↪ Fasten the holding plate with two fixing screws (1) to the printer, see Fig. 5.

↪ Fit the printer with the holding plate onto the corner piece (narrow).

Printer cable fitting

There are 4 cables for connecting the printer. The connection plugs are coded such that incorrect connection is not possible.

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NOTE

The wiring harness would be damaged by sharp edges or hot surfaces.

↪ Do not lay cables over sharp edges or hot surfaces.

↪ Use suitable edge protection if necessary.

↪ Lay cables along existing wiring harnesses.

↪ Connect the USB printer connection cable (5) connection plug to the connection socket (6) of the printer, see Fig. 6.

↪ Connect the USB printer connection cable (5) USB plug to the connection socket (4) of the USB extension cable (3), see Fig. 6.

↪ Lay the USB extension cable to the CPU-1 electronic module (8N1) (1), see Fig. 6.

↪ Connect the USB extension cable plug to the connection socket ST15 on the CPU-1 electronic module (8N1), see Fig. 7.

↪ Connect the printer adapter cable (8) connection plug to the connection socket of the printer, see Fig. 6.

↪ Lay the printer adapter cable (8) from the printer to the connector 9X5-3 (9) in the machine, see Fig. 6.

↪ Connect the printer adapter cable (8) connection plug to the connector 9X5-3 (9) in the machine, see Fig. 6.

↪ Connect the printer connection cable (10) connection plug to the connector 9X5-3 (9) in the machine, see Fig. 6.

↪ Lay the printer connection cable from the connector 9X5-3 (9) to the connector 16X5-3 (11) in the machine, see Fig. 6.

↪ Connect the printer connection cable (10) connection plug to the connector 16X5-3 (11) in the machine, see Fig. 6.

Printer activation

- ✍ Log in with a service technician with administrator rights.
- ✍ Select the System menu.
- ✍ Select the Printer submenu.
- ✍ Set the printer parameters.

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el**en-US****Printer (PR)**

Reason: This conversion kit enables the internal printer (PR) option to be installed in models PLW 8615, PLW 8616, PLW 8617, PWD 8626 or PWD 8628.

Parts required

Quantity	Mat. no.	Designation
1	10931290	Printer conversion kit

Parts included

Quantity	Mat. no.	Designation
1	10490720	Printer, PG 86xx

Quantity	Mat. no.	Designation
1	10842950	Mounting bracket
2	01023022	Galvanized raised-head self-tapping screw, 2.9 x 9.5
10	00039000	Cable ties, 99 x 2.5
1	10706650	Printer seal
1	10764210	Paper roll, 58 mm
1	10831140	Printer connection cable 8615/8616/8626
1	10831120	Printer connection cable 8617/8628
1	10721040	Printer adapter cable 8615-8628
1	10953760	USB 2.0 extension cable type A/A
1	10953770	USB 2.0 extension cable type A/A

The TSD will be needed for all service work

The warning and safety instructions in the model-specific TSD must be observed.

Service work must be carried out following the instructions in the model-specific TSD.

Printer (PR) (12N1)

Legend for Fig. 1:

1. PIN 1 position plug ST2 USB 2.0
2. PIN 1 position plug ST1 power supply 5VDC/< 3A
3. Holding clips
4. Thread for retaining screws
5. LED - Operating states
6. Paper feed button
7. Confirmation button for continuing with print after paper roll replacement

Printer technology: Thermal, direct

Operational state	Description/State
Constantly off	Power off
1x flash - Pause ¹⁾	Power on/Not ready
Constantly on	Ready for operation
2x flash - Pause ¹⁾	Fault: Printer head
3x flash - Pause ¹⁾	Fault: Paper end or paper feed fault
4x flash - Pause ¹⁾	Fault: High voltage, low voltage or temperature fault

LED - Operating states

¹⁾ = Pause between the flashing code = 1 s (flashing rhythm = 128 ms on 128 ms off (8Hz))

Parameters	Units	Value
Temperature during operation	°C	-5 to 50
Moisture during operation	% uncondensed relative humidity	20 - 85
Storage/Transportation temperature	°C	-40 to 80
Moisture during storage/transport	% uncondensed relative humidity	10 - 90

Ambient parameters

Component	Material
Cover	Sabic Lexan 243R (GE)
Base	Sabic Cycloy C2950

Material**Printer installation**

Depending on model (900 mm or 1150 mm width), only one of the following procedures can be carried out:

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Continue with **Preparing the printer for installation**.

Wide corner piece opening on unloading side

- ↺ Open the top box paneling service flap (A); see Fig. 2.
- ↺ Release the locks on the bottom service panel (B); see Fig. 2.
- ↺ Remove the bottom service panel (C); see Fig. 2.
- ↺ Remove the top retaining screw (D) and the bottom retaining screw (E) from the wide corner piece; see Fig. 2.
- ↺ Slide out the wide corner piece partially from the machine (F); see Fig. 2.
- ↺ Open the wide corner piece (G); see Fig. 2.
- ↺ Remove the printer opening cover plate on the corner piece.
- ↺ Continue with **Preparing the printer for installation**.

Narrow corner piece removal on unloading side

- ↺ Open the top box paneling service flap (A); see Fig. 3.
- ↺ Release the locks on the bottom service panel (B); see Fig. 3.
- ↺ Remove the bottom service panel (C); see Fig. 3.
- ↺ Remove the retaining screw for the narrow corner piece and fascia (D); see Fig. 3.
- ↺ Lift the fascia slightly and use a piece of wire or a cable tie inserted through the screw hole to secure it to the top box paneling holder.
- ↺ Lift the narrow corner piece slightly (E); see Fig. 3.
- ↺ Remove the narrow corner piece (F); see Fig. 3.
- ↺ Remove the printer opening cover plate on the corner piece.

↪ Continue with **Preparing the printer for installation**.

Preparing the printer for installation

↪ Remove the holding clamps (4) from the printer (3); see Fig. 4.

↪ Install the printer (3) with its seal (1) in the cutout in the wide corner piece (2); see Fig. 4.

Installing the printer in the wide corner piece

↪ Reinstall the holding clamps (4) in the retainers on the printer (3); see Fig. 4.

↪ Press in the holding clamps (4) until they engage in the wide corner piece (2) and thus secure the printer (3) to the wide corner piece; see Fig. 4.

Installing the printer in the narrow corner piece

The printer holding clamps are only required if the printer is installed in the wide corner piece.

For installing the printer in the narrow corner piece, a special mounting bracket is supplied with the kit.

↪ Install the mounting bracket over the printer so that the openings for the connection plug are on the underside.

↪ Fasten the mounting bracket to the printer with two retaining screws (1); see Fig. 5.

↪ Install the printer with the mounting bracket on the narrow corner piece.

Printer cable installation

There are 4 cables for connecting the printer. The connection plugs are coded so that incorrect connection is not possible.

- Printer connection cable.
- Printer adapter cable.
- USB printer connection cable.
- USB extension cable.

NOTE

The wiring harness can be damaged by sharp edges or hot surfaces.

- ↪ Do not lay cables over sharp edges or hot surfaces.
- ↪ Use suitable edge protection as necessary.
- ↪ Lay cables along existing wiring harnesses.

↪ Connect the USB printer connection cable (5) connection plug to the connection socket (6) of the printer; see Fig. 6.

↪ Connect the USB printer connection cable (5) USB plug to the connection socket (4) of the USB extension cable (3); see Fig. 6.

↪ Lay the USB extension cable to the CPU-1 electronic module (8N1) (1); see Fig. 6.

↪ Connect the USB extension cable plug to the connection socket ST15 on the CPU-1 electronic module (8N1); see Fig. 7.

↪ Connect the printer adapter cable (8) connection plug to the connection socket of the printer; see Fig. 6.

↪ Lay the printer adapter cable (8) from the printer to connector 9X5-3 (9) in the machine; see Fig. 6.

↪ Connect the printer adapter cable (8) connection plug to connector 9X5-3 (9) in the machine; see Fig. 6.

↪ Connect the printer connection cable (10) connection plug to connector 9X5-3 (9) in the machine; see Fig. 6.

↪ Lay the printer connection cable from connector 9X5-3 (9) to connector 16X5-3 (11) in the machine; see Fig. 6.

↪ Connect the printer connection cable (10) connection plug to connector 16X5-3 (11) in the machine; see Fig. 6.

Printer activation

- ↗ Log in as a service technician with administrator rights.
- ↗ Select the System menu.
- ↗ Select the Printer submenu.
- ↗ Set the printer parameters.

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