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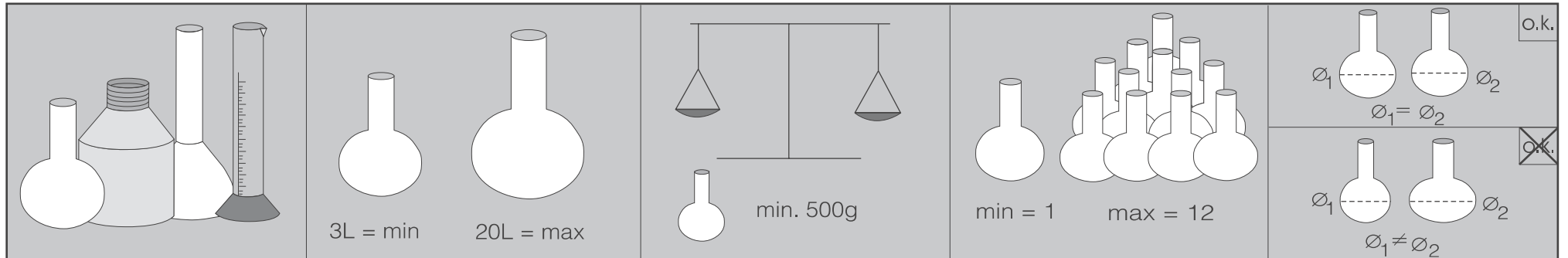
⚠ Please observe the warning and safety instructions given in the operating instruction manual supplied with your washer disinfector, as well as your laboratory glassware manufacturer's instructions.
 Glass broken during the cleaning process can cause serious injury. Cleaning results should be tested properly. Do not rely on visual checks only.

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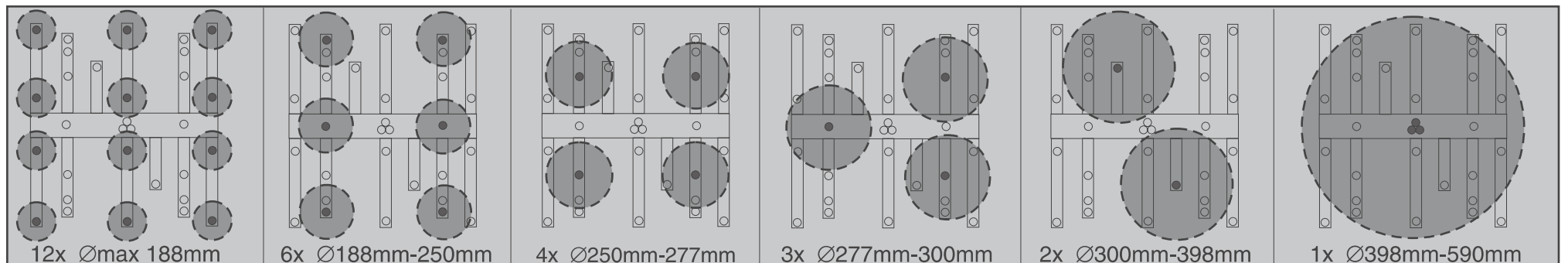
The E 957 should only be used for processing laboratory glassware which has been declared as suitable for machine processing by the glassware manufacturer.








Laboratory glassware, which because of its design cannot be flushed through properly, will not be sufficiently cleaned and disinfected in this mobile unit.

This mobile unit is not designed to be continually rearranged for different loads.

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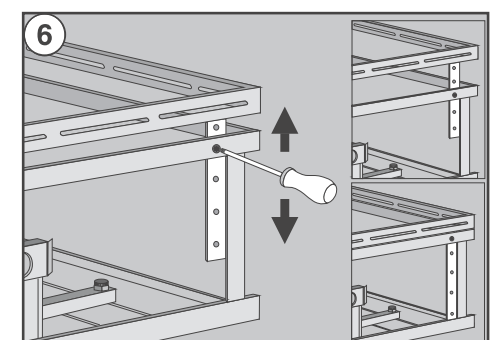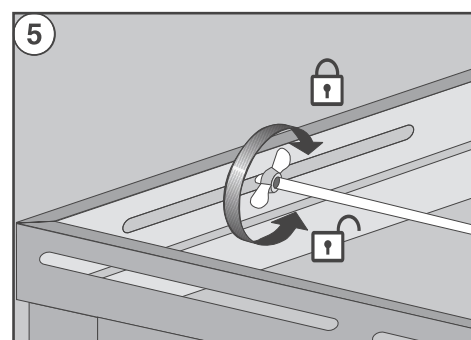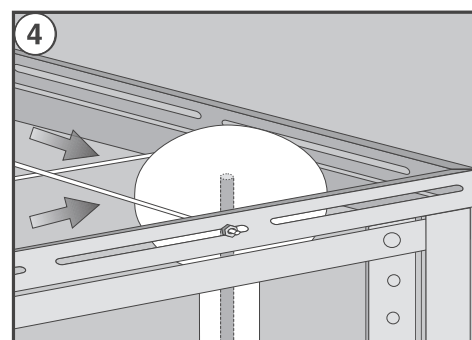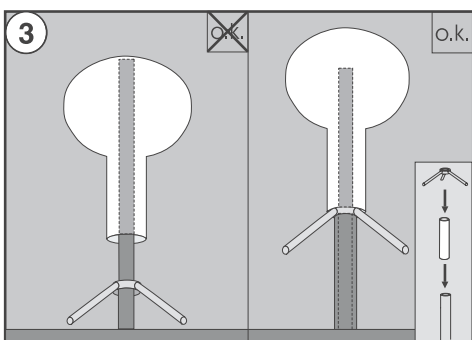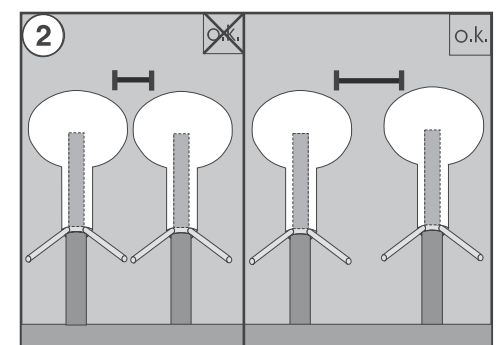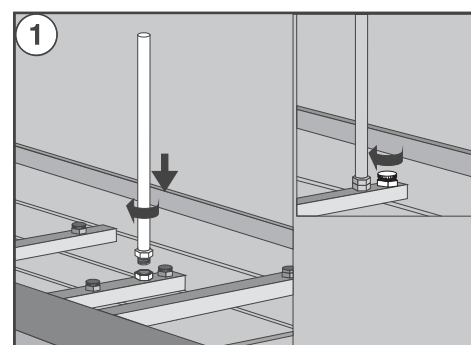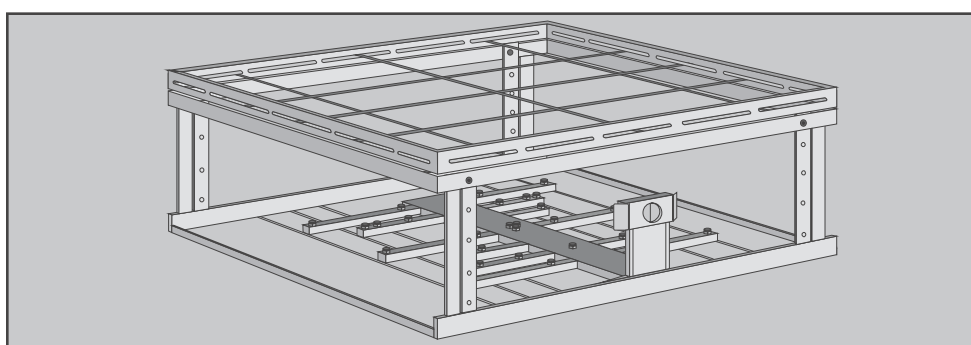Handhabungshinweise / Setting up the mobile unit

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- Screw the jets into the positions required. Unused positions should be plugged using blanking screws to ensure sufficient water pressure is maintained (1).
- Place the laboratory glassware over the jets. Make sure there is sufficient space between each item of glassware (2).
- Spacers can be used to hold the bottom of flasks etc. off the top of the jets (3).
- Secure the laboratory glassware in position with rods (4).
- Tighten the wing nuts to secure the rods and hold in place with hex. nuts (5).
- If the laboratory glassware is too tall to secure at its widest diameter you will need to use spacers or adjust the height of the mobile unit's frame (6).

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